ACM

USER'S MANUAL MANUAL DE USUARIO MANUEL DE L'UTILISATEUR BEDIENUNGSANLEITUNG

EN - ES - FR -DE

PRECISION SCALES

BALANÇA PRECISIÓN

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V.1.1
20180602

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EN / CONTENTS

Warming and safety Using	
I	First part: Summarize 5
II	Second part: Installation 7
III	Third part: Basic operation 8
	BASIC WEIGHING FUNCTION 9
	CALIBRATION / ADJUSTMEN 9
IV	Fourth part: Operate Application 9
	COUNTING FUNCTION 9
	PERCENTAGE WEIGHING 9
	ACCUMULATING 9
V	Fifth part: Basic Function Setting 9
	BACKLIGHT SETTING 9
	POWER ON SELF TEST SETTING 9
	BUZZER SETTING 9
	TURN ON/OFF UNITS 9
VI	Sixth part: Machine Weighing Configuration Setting. 11
VII	Seventh part: restore factory setting 12
VIII	Eighth part: machine detail specification 12
IX	Ninth part: proper care and maintenance 13

We ware try the best to ensure the veracity of operationg manual, but we didn't take responsibility for printing or description mistake.

We has right update the machine looking and performance without noticing the costumer.

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WARNING AND SAFETY USING

SAFETY

For avoiding damage, please read all operating instructions carefully before use.

- ⚠ Don't use your machine under dangerous working circumstance
- ⚠ Cut off the power if machine will turn off for more than one week
- ⚠ Turn off the machine and cut off the power before or after connection with other equipments.
- ⚠ Strong magnetic field and static electricity can have an adverse effect on weighing sensor. When disturbance disappears, the machine will work well again.

WARNING

- » All our parts is the most suitable parts for machine.
All modification or using unauthorized parts for machine need to be confirmed before using.
All modification needs to be taken responsibly.
- » Do not open the machine housing. Machine will not have guarantee service if security label broken.

FIRST PARTH: SUMMARIZE

1. UNPACKING

After unpacking the machine, please check machine has any visible damage.

Please keep the original box and packing material for storing material for storing machine when not in use or send back for repairing.

Before packing the balance, please cut off all power and cable.

2. PACKING LIST

Main machine	Adapter	Pan	Operating Manual	Calibration mass
1pc	1pc	1pc	1pc	1pc

3. INSTALLATION

When select the location for install machine, please keep these tips in mind:

- Do not put machine close to central heating or sunshine and airflow way. (Opening door window).
- Do not exposure machine to extreme heat or cold. Keep scale in a clean, dry location. Dust, dirt and moisture can accumulate on the weighing sensor.
- Install machine on a flat and level surface, free from vibration and drafts, free from corrosive and strong magnetic field, as they can have an adverse effect on the weighing sensors.

4. WARM UP FOR MACHINE ADAPT TEMPERATURE

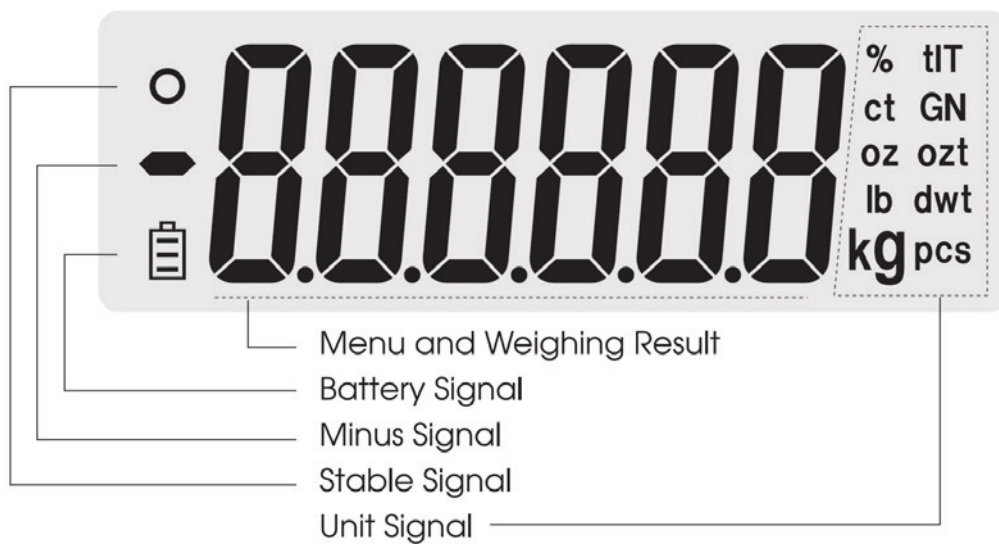
When move machine from high temperature place to low temperature place (or inversely), please keep machine final place for two hours and then turn on to warm up (warm up time refer to the specification list), as the machine will proportion the room temperature.

5. KEY EXPLANATION

Key symbol	
U	<p>UNIT KEY</p> <p>A: Select Unit. B: Status 1: Enter into menu with (TARE) KEY C: Status 2: Setting Key D: Status 3: Shortcut key of accumulating function. Hold and press this key to enter into accumulating function.</p>
	<p>CAL KEY</p> <p>A: When normal weighing, short press CAL KEY will zeroing. B: Press and hold CAL KEY for 5 seconds will enter into CALIBRATION. C: When setting the parameter, press CAL KEY can minus one value. D: CONFIRM KEY.</p>
	<p>MODE KEY</p> <p>A: Short press MODE KEY will enter into Counting Function. B: Move the flash on digit to left.</p>
	<p>TARE KEY</p> <p>A: Tare the weight B: Move the flash on digit to left.</p>
	<p>Press and Hold TARE and UNIT KEY at the same time for 5 seconds enter into setting menu.</p>

NOTE: The buzzer sound different when long press or short press the key.

6. DISPLAY LCD

II SECOND PART: INSTALLATION

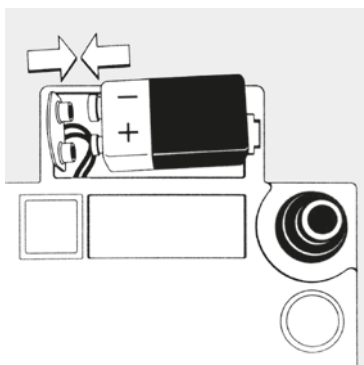
The machine with round weighing pan

- Put weighing pan on the pillar which is in the middle of machine

The machine with square weighing pan

- Put the weighing pan on bracket.

Use dry battery / Rechargeable battery (Optional)

The dry battery or rechargeable battery is not on machine's packing list.

- ⚠ Only normal or universal 9V dry battery will be available for machine.
- Lie down the machine at side.
- Open the battery cover box cover.
- Connect and put 9V dry battery.
- Confirm the positive and negative correctly.
- Close the battery box: screw the battery box cover down to the machine.

⚠ The used battery is recycled. According to the waste disposal law.

Adjust machine Level

The machine needs to be adjusted every time when the installation location is changed. Moving the two back screws slowly to adjust the level.

- Counterclockwise rotate the two back screws to the right position.
- Rotate the screws as shown in the photo until the bubble is in the middle of the level device.

- Clockwise rotate the two back screws until they touch the support.

> Under normal circumstances, adjusting the level may need several times to reach the suitable position.

III THIRD PART: BASIC OPERATION

1 BASIC WEIGHING FUNCTION

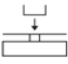
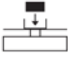
PREPARATION

Turn on machine: Press (ON/OFF) Key

WARM UP TIME:

For making sure the weighing result correct, different type machine need different warm up time to reach the required operating temperature. please refer the specifications list to know the correct warm up time.

INSTANCE:

Basic Weighing (The machine warm up)		
Key (Order)	Step Explanation	LCD Screen Display
	1. Zero Stable	0.00g
	2. Put on container on weighing pan (Example: 100g)	100.00g
	3. Press Tare key for tare the container weight	0.00g
	4. Put sample in container (Example: 200g)	200.00g

2 CALIBRATION / ADJUSTMNET

Request

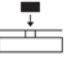
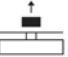
The machine only can be calibrated/Adjusted in the following cases:

- Nothing on weghing pan, no tare, weighing signal stable.

Machine will have ERROR SIGNAL if does not have above premise.

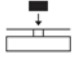
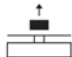
If all premise available, the machine will display required will display required weight mass value for calibration.

INSTANCE: (Example: A200)

(1) External Span Calibration Instance		
Key (Order)	Step Explanation	LCD Screen Display
Press and hold  key	1. Scale will ask for the calibration mass.	200.00g
	2. Put a calibration mass on weighing pan (Example: 200g)	200.00g
	3. Wait untill the indication stops flashing.	0.00g

(2) External Linearity Calibration Instance

(Please DO NOT linearity calibrate the machine if you don't have match weight mass)

Key (Order)	Step Explanation	LCD Screen Display
Short press  key	1. Machine Tare	0.00g
Press and Hold  key for 5 seconds.	2. Display Span CAL signal Flash 200g signal	200.00g
Release  key, Press and Hold  key for 5 seconds.	3. Display Linearity CAL signal Flash 100g signal	100.00g
	4. Put required weight mass on pan Display 100g after 5 seconds.	100.00g
	5. Take weight mass away (Linearity calibration finished)	0.00g

Note: The grey color words explain the signal's meaning which flash on the window.

IV FOURTH PART: OPERATE APPLICATION

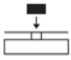
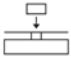
APPLICATION SETTING

COUNTING

PURPOSE: Use this function can calculate the quantity with total weight divide by signal weight.

INSTANCE: (With know the sample's quantity but unknown the unit weight)

Basic Weighing (The machine warm up)		
Key (Order)	Step Explanation	LCD Screen Display
Short Press  key	1. Enter into counting function and display flash the number of quantity. Short press  key to cycle the sample quantity setting: 10, 20, 50, 100, 150, 200, 250, 500, 1000pcs.	
	2. Put any unknown numbers of pieces on pan. (Example: set 20pcs)	20 pcs
Short Press  Key	3. The machine start to calculate and display quantity: 20pcs	20 pcs
	4. Take sample away from pan	0 pcs
	5. Put any unknown numbers of pieces on pan and will display a count. (Example: 100pcs)	100 pcs
Short Press  Key	6. Exit the counting function	0.00 g

Basic Weighing (The machine warm up)		
Key (Order)	Step Explanation	LCD Screen Display
	1. Zero Stable	90.00g
Short Press U Key	2. Put container on weighing pan (Example: 100g)	100.00%
	3. Take sample away and put any other sample on pan. Remove the reference sample and add the unknown sample to determine its relative weight and percentage.	79.00%
Short press [UNIT] key	4. Exit the percentage Measurement	0.00g

V FIFTH PART

PURPOSE

Operation can set machine basic function by selecting parameter in menu.

POWER ON SELF TEST SETTING (POST)

User can enter into the menu to turn on or off the Power On Self Test of voltage and temperature.

BUZZER ON/OFF SETTING

Operator can turn on/off the buzzer sound by setting menu.

BASIC FUNCTION SETTING

LCD Basic setting Instance

Key (Order)	Step Explanation	LCD Screen Display
Press and Hold [TARE]&[UNIT] Key for 5 seconds at the same time.	1. When display the menu, enter cod:101 The code of basic setting menu is C00101. The enter way is: Press [MODE] key to move the digit, press the [TARE] key to circle and set the number.	C00 10 1
Short Press [UNIT] Key	2. Enter into back light setting menu, display BL and flash ON (turn on) Press the [TARE] key or [CAL] key can circle and select the backlight: ON / turn on), OFF (turn off), AUT(turn on or off the backlight automatically).	BL - ON
Short Press [UNIT] key.	3. Enter into Power On Self Test menu, display POST and flash O (0 means not show the test). Short press [TARE] key can select 0 (not show) or 1 (show the test).	POS
Short Press [UNIT] key.	3. Enter into Buzzer setting menu, display BEP and flash ON (turn on) Short press [TARE] key can select ON (turn on) or OFF (turn off) the buzzer.	100.00g
Short Press [UNIT] key.	4. Set the machine successful and return to the main menu.	C00000
Short Press [UNIT] key.	5. Return to standby	

The green color parts is the following operation after Step 1-2 backlight setting.

NOTE: The grey color words explain the signal's meaning which flash on the window.

Turn On/Off the units

Operator can turn on or off the unit to display or hide the relative weighing units.

INSTANCE

Key (Order)	Step Explanation	LCD Screen Display
Press and Hold [TARE]&[UNIT] Key for 5 seconds at the same time.	1. When display the menu, enter cod:106 The code of basic setting menu is C00101. The enter way is: Press [MODE] key to move the digit, press the [TARE] key to circle and set the number.	<i>C00 106</i>
Short Press [UNIT] Key	2. Enter into the menu, display unit: ct and flash ON (turn on)	<i>ct ON</i>
Short Press [TARE] key.	3. Display unit: ct and FLASH (turn off) (Set OFF/ON with short press repeatedly)	<i>ct OFF</i>
Short Press [MODE] key.	4. Enter into Buzzer setting menu, display BEP and flash ON (turn on) Short press [TARE] key can select ON (turn on) or OFF (turn off) the buzzer.	<i>o2 ON</i>
Short Press [TARE] key.	5. Display unit: oz and flash OFF Repeat the step 4-5 can circle and set follow units: ct, oz, ozt, lb, dwt, tIT, kg, % The default setting is all units was turn on.	<i>o2OFF</i>
Short Press [UNIT] key.	6. Set the machine successful and return to the main menu.	<i>C00000</i>
Short Press [UNIT] key.	7. Return to standby	

NOTE: The grey color words explain the signal's meaning wich flash on the window.

VI SIXTH PART

Machine Weighing Configuration Setting

Purpose

Operator can set the machine basic weighing config to change the weighing capability to reach different required.

Zeroing Range Setting

Operator can increase or decrease tracking range for they need.

Tracking Range Setting

Operator can increase or decrease tracking range for they need.

Sensitivity Level Setting

Operator can adjust the weighing response time by increase or decrease the level. Level 1 is the slowest weighing resopnse speed and level 3 is the fastest (Default and recommend setting: Level 2)

Speed Level Setting

Operator can adjust the weighing response time by increase or decrease the level. Level 1 is the slowest weighing response speed and level 3 is the fastest (Default and recommend setting: Level 2)

Anti-Vibration level Setting

Operator can adjust the weighing response time and anti-vibration strength by oncrease or decrease the level.

The higher level comes with higher anti-vibration. Level 1 has fast weighing speed and weak anti vibration. Level 7 has strong anti vibration and low weighing speed.

INSTANCE		
Key (Order)	Step Explanation	LCD Screen Display
Press and Hold [TARE]&[UNIT] Key for 5 seconds at the same time.	1. Display the menu and enter into the code: "100". The weighing config: C00100, enter way: Press [MODE] key to move the digit, press [TARE] key to circle and set the number.	<i>C00 100</i>
Short Press [UNIT] Key	2. Enter into the zeroing range setting. Short press [TARE] key or [CAL] key can select zeroing range from 0.0 to 6.0	<i>2Er00.0</i>
Short Press [UNIT] key.	3. Enter into tracking range setting. Short press [TARE] key or [CAL] key can select tracking range from 0.0 to 6.0	<i>SLD405</i>
Short Press [UNIT] key.	3. Enter into Sensivity Level Setting Short press [TARE] key or [CAL] key can select Sensitivity Level from 1 to 6. Totally six levels.	<i>SENS-3</i>
Short Press [UNIT] key.	3. Enter into Speed Level Setting Short press [TARE] key or [CAL] key can select Speed Level from 1 to 3. Totally three levels.	<i>SPEED2</i>
Short Press [UNIT] key.	3. Enter into Anti-vibration Level Setting Short press [TARE] key or [CAL] key can select anti-vibration Level from 1 to 7. Totally seven levels.	<i>FILL-4</i>
Short Press [UNIT] key.	4. Setting successful and return to the main menu.	<i>C00000</i>
Short Press [UNIT] key.	5. Return to standby	

The grew color parts is the following operation after Step 1-2 Zeroing Setting.

NOTE: The grey color words explain the signal's meaning wich flash on the window.

VII SEVENTH PART: UNIT SWITCHING

Press [UNIT] key, the weighing unit will cycle between the different weighing units with each press of the button. The balance will default the last unit used when turned on the next time.

LED/HD/DJ Series

Unit Signal	Unit	Unit Exchange Rate
g	Gram	1
ct	Carat	5
oz	Ounce	0.03527396200
ozt	Troy Ounce	0.03215074700
dwt	Pennyweight	0.64301493100
GN	Grains	15.43235835000
lb	Pound	0.00220462260
N	Newton	0.00980654189
tlT	Taiwan Tael	0.02666666000
kg	Kilogram	0.00100000000

VIII EIGHTH PART: MACHINE DETAIL SPECIFICATION

Device	Weighing Range (g)	Readability (g)	Repeatability (g)	Linearity (g)	Operate Temp (°C)	Pan Size (mm)	Housing Size (LxWxH)(mm)	Warm-up Time (min)
ACM 2	200	0.01	±0.01	±0.02	10–35	Ø 128	260X175X60	10–20
ACM 3	300	0.01	±0.01	±0.02	10–35	Ø 128	260X175X60	10–20
ACM 5	500	0.01	±0.01	±0.02	10–35	Ø 128	260X175X60	10–20
ACM 6	600	0.01	±0.01	±0.02	10–35	Ø 128	260X175X60	10–20
ACM 10	1000	0.01	±0.01	±0.02	10–35	Ø 128	260X175X60	10–20
ACM 20	2000	0.01	±0.02	±0.03	10–35	Ø 128	260X175X60	10–20
ACM 30	3000	0.01	±0.01	±0.02	10–35	145X160	145X160	10–20
ACM 50	5000	0.01	±0.01	±0.02	10–35	145X160	145X160	10–20
ACM 60	6000	0.01	±0.01	±0.02	10–35	145X160	145X160	10–20

IX NINTH PART: PROPER CARE AND MAINTENANCE

REPAIR

Only trained technician was authorized to repair the problem machine.

CLEAN

- Pull out the adapter from electrical outlet and cable from machine.
- Use soft cloth with neutral cleanser to clean the machine housing.
- Dry the housing with soft cloth and then take out the weighing pan and wash it.
- When take up the weighing pan and bracket, make sure that don't broken the weighing system.

⚠ Do not let the liquid flow into machine.

⚠ Do not use the caustic cleanser.

WASH STAINLESS STEEL SURFACE

Use soft cloth or sponge to clean all stainless steel parts need to clean often and completely. Only home appliances cleanser available for clean the stainless parts.

Wipe up the stainless steel parts surface first, wash up all leftover second and then dry it. Oil the stainless steel surface if necessary.

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marca propiedad de | est une marque de | trade mark propriety of | die Marke ist Eigentum von:
marca propiedad de | est une marque de | trade mark propriety of: